METTLER TOLEDO

InPro 3100 (i) pH Electrode

Instruction Manual

cs	Návod k obsluze	2
da	Brugsanvisning	6
de	Betriebsanleitung	10
en	Instruction Manual	14
es	Manual de instrucciones	18
fr	Instructions d'utilisation	22
fi	Käyttöohje	26
hu	Használati utasítás	30
it	Istruzioni per l'uso	34
ja	取扱説明書	38
ko	사용 매뉴얼	42
nl	Gebruikershandleiding	46
pl	Instrukcja obsługi	50
pt	Manual de instruções	54
ru	Инструкция по эксплуатации	58
sv	Bruksanvisning	62
th	คู่มือคำแนะนำ	66
zh	说明书	70

InPro 3100 (i) pH elektroda

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Obsah

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
Instruction Manual

Contents

1	Introduction	15
2	Safety instructions	15
3	Product description.....	15
4	Installation and commissioning	16
5	Operation.....	16
5.1	Calibration of the electrode and pH transmitter	16
5.2	Sterilization of electrode	16
5.3	Temperature sensor of InPro 3100 (analog version).....	16
6	Maintenance	17
7	Trouble-shooting (sluggish response time, insufficient sensitivity etc.)	17
8	Disposal.....	17
9	Guarantee.....	17

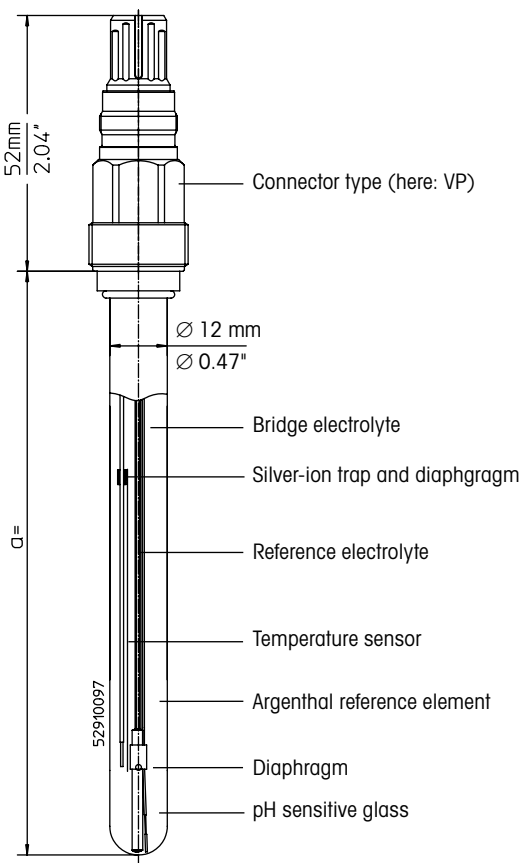InPro and ISM are registered trademarks of METTLER TOLEDO Group in Switzerland, the USA, the European Union and a further five countries.

1 Introduction

The METTLER TOLEDO InPro™ 3100 (i) pH electrodes are low-maintenance, autoclavable and pressure-resistant measuring electrodes with a gel-type reference electrolyte. The reference element incorporates a silver-ion trap to prevent silver sulfide precipitation on the diaphragm. In case of InPro 3100, a temperature sensor is integrated, mainly for automatic temperature compensation possibilities of the pH signal during calibration and operation.

The InPro 3100 (i) electrodes are mainly designed for the measurement of pH value in biotechnical processes at moderate pressures (<6 bar). It can be used under industrial fermentation, pilot and laboratory conditions, and is particularly intended for those processes where intermediate sterilization or autoclaving is indispensable.

Please read through these operating instructions carefully before commissioning, in order to ensure trouble-free use.

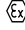
2 Safety instructions

The InPro 3100 (i) electrodes are intended solely for the purpose described in §1 above. The electrode consists partly of glass and should be handled carefully. If the electrode is to be cleaned or calibrated using acid or alkaline solutions, safety goggles and gloves should be worn. Local safety regulations are applicable to the handling of the electrode. Please follow the guidelines enclosed for an installation in the Ex-zone.

We recommend that you only operate the electrode in combination with original parts from METTLER TOLEDO. Operation and maintenance should be carried out only by trained personnel and staff who have read and understood the operating instructions.

3 Product description

The inscription on each electrode contains the following information:

METTLER TOLEDO	Manufacturer of the electrode
InPro™ 310XiUD/SG/yyy/zzz	Type designation; X = membrane glass; i = ISM Digital; UD = upside down/SG = ancillary electrode/yyy = rod length in mm/zzz = type of temperature sensor
combination pH	Type of electrode
pH 0 ... 14	pH measurement range
0 ... 80 °C / 140 °C (130 °C)	Temperature range for operation/sterilization (for InPro 3100UD)
Order No. 52 00X XXX	Order number
 SEV14 ATEX 0168 X	Identification and number of certificate
 IECEx SEV 14.0025X	Identification and number of certificate
 CML 22 UKEX 2108X	Identification and number of certificate
IS/I, II, III/1 / ABCDEFG/T6	FM approved
See instruction manual!	Instruction (Observe the operating instructions)

In addition each electrode is allocated a serial number on the plug-in head to enable identification and traceability.

4 Installation and commissioning

1. On unpacking, check the electrode for mechanical damage of the pH sensitive glass and shaft. Report any damage immediately to your METTLER TOLEDO supplier.
2. Remove the watering cap and briefly rinse the electrode with deionized water. After rinsing, the electrode should only be dabbed dry with a tissue. Do not rub the pH sensitive glass, since this can lead to electrostatic charging and sluggish response times.
3. Check the space behind the pH sensitive glass for the presence of any air bubbles and remove same by gently swinging the electrode in a vertical plane.
4. Please refer to the instruction manual of the relative housing in order to install the electrode.
5. Connect the electrode to the pH transmitter, using a cable with the appropriate connector (VP connector for InPro 3100). Please refer to the wiring and terminal diagram accompanying the cable.
6. Intelligent ISM[®] electrodes such as InPro 310Xi enable "Plug and Measure" and extended diagnostics. For installation, commissioning and operation of the ISM system please observe the instructions for the transmitter, module and cable.

5 Operation

5.1 Calibration of the electrode and pH transmitter

For the InPro 3100(i) electrode a 2-point pH calibration is recommended. Before calibration, first remove the watering cap. To speed up the response time of the electrode after long transport and storage periods, and before calibration, dip the electrode into buffer pH 7, buffer pH 4 and buffer pH 7 again, each time for about one minute. The electrode is then dipped successively into two different buffer solutions with given pH reference values and the transmitter calibrated to these buffer values. Normally you start with buffer pH 7 for determining the 0-point and then buffer pH 4 or 9 for determining the slope. In the case of microprocessor-based transmitters with buffer recognition, the order of the two chosen buffer solutions is irrelevant. Whenever possible, a temperature sensor (for InPro 3100 the integrated one) should be connected to the transmitter and "automatic temperature measurement" selected there. Please refer to the instruction manual of the transmitter for further details.

5.2 Sterilization of electrode

The InPro 3100(i) electrode can be sterilized either through autoclaving (remove red protection cap) or through in-situ sterilization with steam or superheated fermentation medium. Sterilization temperatures above 140 °C are not allowed, and excessive sterilization time (for instance $t > 30$ min at 140 °C) could notably shorten the life time of the electrode. Repeated sterilization and/or autoclaving of the electrode can lead to a slight coloration of the electrolyte. This color change does not effect the functioning of the electrode.

5.3 Temperature sensor of InPro 3100 (analog version)

The integrated Pt100/1000 temperature sensor in the InPro 3100 is primarily intended for automatic temperature compensation of the pH signal and **not** for any accurate and safe temperature indication or control of the process temperature. If you anyhow want to try this, check that your transmitter has the required temperature indication accuracy and a free outlet for the signal. Furthermore, a calibration of the temperature signal might be necessary.

6 Maintenance

1. Carefully rinse the electrode tip and diaphragm with deionized water after each production cycle. Under no circumstances must measuring solution be allowed to dry on these parts!
2. When the electrode is not in operation, store it with electrode tip and diaphragm well submerged into 9823/3M KCl or 9816/Viscolyt™ electrolyte.
3. If the electrode is stored mounted in its housing, the same rule as in 2) above applies, but the storage electrolyte has to be slightly modified, with 2 parts of buffer solution pH 9.2 to 10 parts of electrolyte, this in order to avoid corrosion of housing parts.
4. If an electrode is stored dry for a few days by mistake, let it soak in the normal storage electrolyte for several hours before use.
5. Occasionally check the connector for possible traces of moisture. If necessary, clean well with deionized water or alcohol, and afterwards dry carefully.

7 Trouble-shooting (sluggish response time, insufficient sensitivity etc.)

The glass membrane can be reactivated using a HF solution (order no. 51 319 053). The electrode tip (only pH sensitive glass part) is immersed into the HF solution for a period of about 5 minutes, is then immediately and thoroughly rinsed with deionized water and thereafter stored in buffer solution pH 4 (see point 6 above) up to the diaphragm for 12 hours. Possible protein precipitation and blockage of the diaphragm can be treated with our special electrode cleaner (order no. 51 340 068). However, use of the diaphragm cleaner (order no. 51 340 070) is normally not necessary for this electrode type, as the built-in silver-ion trap prevents the formation of silver sulfide contamination.

Also, check the connector for possible traces of moisture. If necessary, clean well using deionized water or alcohol, and afterwards dry carefully.

8 Environmental protection

Waste electrical products should not be disposed of with household waste. Please recycle where facilities exist. Check with your Local Authority or retailer for recycling advice.

9 Guarantee

On manufacturing defects, 12 months after delivery.

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 本表依据SJ/T 11364的规定编制。本产品符合以下标志规范：
 Table composed in accordance with SJ/T 11364 (CN). This product is bearing the following symbol:

○ : 表示该有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下
 ○ : Indicates that the content of the hazardous substance in all homogeneous materials of the part is below the limit specified in GB / T 26572

x : 表示该有害物质至少在该部件的某一均质材料中的含量超出GB/T 26572规定的限量要求。
 x : Indicates that the content of the toxic substance in at least one of the homogeneous materials of the part exceeds the limits specified in GB/T 26572.

产品中有毒物质的名称及含量 Toxic and hazardous substance name and containment in product						
部件名称 Part Name	有毒有害物质或元素 Toxic and hazardous substances					
	铅 Lead (Pb)	汞 Mercury (Hg)	镉 Cadmium (Cd)	六价铬 Hexavalent chromium (Cr6+)	多溴联苯 Polybrominated biphenyls (PBB)	多溴二苯醚 Polybrominated diphenyl ethers (PBDE)
插头 Elect. Connector	x	○	○	○	○	○
电路板 PCBA	x	○	○	○	○	○

本表依据SJ/T 11364的规定编制。本产品符合以下标志规范。
Table composed in accordance with SJ/T 11364 (CN). This product is bearing the following symbol:

○：表示该有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下
○：Indicates that the content of the hazardous substance in all homogeneous materials of the part is below the limit specified in GB / T 26572

x：表示该有害物质至少在该部件的某一均质材料中的含量超出GB/T 26572规定的限量要求。
x：Indicates that the content of the toxic substance in at least one of the homogeneous materials of the part exceeds the limits specified in GB/T 26572.

For addresses of METTLER TOLEDO
Market Organizations please go to:
www.mt.com/contacts

METTLER TOLEDO Group
Process Analytics
Local contact: www.mt.com/contacts

Subject to technical changes
© 08/2022 METTLER TOLEDO
All rights reserved
Printed in Switzerland. 52 002 293 C

Management System
certified according to
ISO 9001 / ISO 14001

**UK
CA**₂₅₀₃

CE₁₂₅₈

EAC

www.mt.com/pro

* 5 2 0 0 2 2 9 3 C *